

MesoSPIM Part 2: Image acquisition

This guide explains the steps required to acquire a z-stack on the MesoSPIM lightsheet microscope.

E mesoSPIM Main Wind File View Movement ETL Scripting Log Parameters Alignment Sample Translation Focus Increment Increment LOAD UNLOAD 500 µm 500 µm Rotation STOP Increment MOVEMENT 0.00° X 1182.9 µm Y 5499.3 µm Z 5593.7 µm Angle 0.0° Focus 3000.0 µm Controls 405-488-561-640-Quadrupleblock Right 10 % 488 nm Run elected Acquisitior Run Acquisition List Live 0% Adjust levels O av: none Total Progress 0% 4, 1024) Z_start Planes Rot F_start Laser Intensity Filter Zoom hutterconfi Folder Filename X_pos F end Empty-Al... 1x 100 10 10 Stack 0 0 488 nm Left tmp one.raw Add Row Duplicate Row Delete Row Move Row Up Save Table Tiling Wizard Image Processing Wiz Mark Current XY Mark Rotation Preview Selected Row Delete all Rows Move Row Down Load Table Filename Wiz Focus Tracking Wiz Mark Focus Mark ETL values Preview: Move to Z start Predicted time (HH:MM:SS): 00:00:02 Set Folders Mark State Mark ALL

Written By: José Maria Mateos Melero

INTRODUCTION

In this guide for the ZMB MesoSPIM lightsheet microscope you learn how to locate and focus a sample as well as setting up and acquiring a z-stack.

For a full documentation about the microscope please, visit the project webpage: http://MesoSPIM.org

Step 1 — Find and center the sample



- Press "live" to get the first image and stop.
- Default settings: Right laser illumination; Zoom: 1; 488 nm; Filter 405-488-561-640-quadrupleblock filter. Change this in the corresponding boxes if you need other settings.
- Click "Adjust Levels" for adjusting the histogram.
- Use the Y arrows to move the sample up or down (Y).
- Use the X arrows to move the sample left or right (X).
- Use the Z arrows to bring the sample into the plane of the lightsheet.
- It is recommended, after some movements, to stop the scan and check the position of the cuvettes to avoid collisions between the inner and outer cuvettes.

Step 2 — Focusing the sample



- Manual Focus: use the Focus arrows to focus the lightsheet. For a fine focus, change the focus "Increment" to 20 um.
- AUTO Focus: click on the "AUTO Focus" to launch an optimizer that lets you automatically search for the best focus within a given range from the current position.
- Click on "RUN"
- Once the autofocus has run a new window will appear showing the optimization results. If you are satisfied with the new focus click "Accept".

Step 3 — Adjustment of ETL parameters

III mesoSPIM Main Window	mesoSPIM Main Window	
File View	File View	
Movement ETL Scripting Log Parameters Alignment	Movement ETL Scripting Log Parameters Alignment	
Tunable lenses Left Right Freeze galvos Offset 2.394 \ 2.492 \ 3.400 \ 2.492 \ 4.400 \ 4.	Tunable lenses Left Right Offset 2.494 \ 2.492 \ AmpLieu 0.898 \ 0.804 \ Increment 0.010 \ Choose ETL config file Save current parameters to csv Config/etl parameters.csv	

- Click "Choose ETL config file", select your ETL parameters file. The file is also specific for the clearing method you are using.
- Manual ETL adjustment with sample : For the ETL optimization press on "AmpL = 0" to set the ETL amplitude to 0.
- Hover with the mouse arrow over the corresponding offset value and use the mouse wheel to adjust it. Find the waist in the camera image and bring it to the center.
- Press again on "AmpL = 0" to set the value back to the stored value.
- Press "save current parameters to csv".
- **ETL parameter optimization**: It best performs using a highly transparent sample with sparse but evenly distributed labeling, such as vasculature.
- Manual method without sample using a parked galvo and scattered laser beam: <u>https://www.youtube.com/watch?v=dcJ9a7VA...</u>

Step 4 — Z-stack acquisition



- In the "mesoSPIM Acquisition Manager", click on "Stack 0", followed by "Mark ALL"
- Saving data: Click on "Folders" and select the Network drive X:\
- Press live and use the Z arrows to select the beginning of the stack.
- Click on the "M" button to the left of the "Z_start" to record this position.
- Use the Z arrows to localize the end of your stack. Then press the "M" button to the left of "Z_end".
- Insert the step size in micrometers. The number of planes will be automatically calculated.
- Click "Filename Wizard". In the appearing menu select "save as" BigDataViewerfile: h5; give a name to the file and press "finish".
- Click "Run Acquisition List"