

Olympus ScanR HCS (Irchel) - Part3: Log off / Shut down

How to log off / shut down the Widefield - Olympus ScanR HCS (Irchel) located at the Center for Microscopy and Image analysis, UZH, Room Y24-F-14.

Written By: Joana Raquel Delgado Martins




Step 1 — Log off / Shut down the computer and microscope board



- Remove and clean your sample.
- Make sure the desk and microscope are as clean and tidy as you would like to find them.
- Close ScanR. Check in the booking system for the next booking.
- If there is another booking shortly after you, log off the PC.
- If you are the last user of the day proceed to the next step.

Step 2 — Shut down

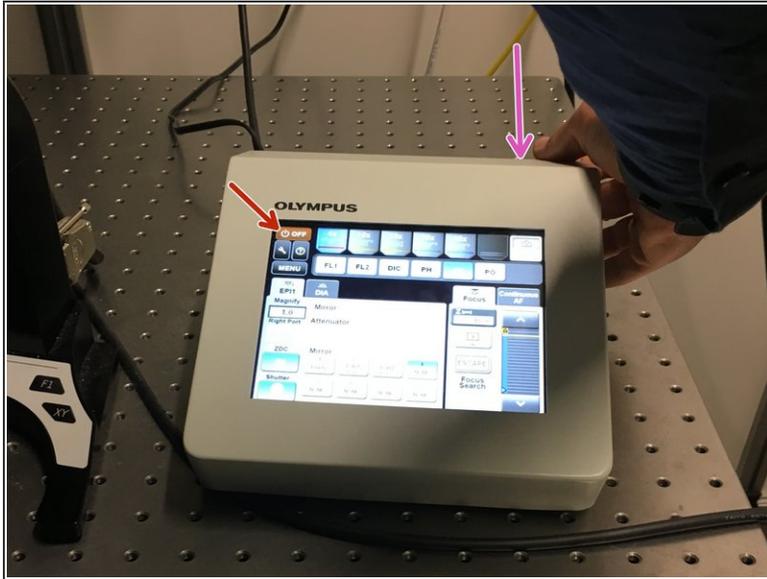

- First switch off the microscope board by clicking "OFF" on the touch screen.
- Switch it off at the back as well.
- Shut down the PC.
- Switch off the **left power** switch on the top shelf.