

# Leica LASX Auto-Save and How to Handle the Resulting Files

This guide shows how to use the auto-save functionality in the Leica LAS X software and how to open the resulting files.

Written By: Dominik Hänni

,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,		Risu — 0				x	
		<ul> <li>Project settings:</li> </ul>			<b>*</b> ()	^	
' I 650		Next project name :	3D_Time_TIF				
		Auto-Save :	ON				
		Base data folder:	\\files.core.uzh.ch\zmbstaff-equipment\t	ransfer\DH_JM_Autos	Browse		
		Current working folder:			Browse		Sec. 1
		Current data path :	\\files.core.uzh.ch\zmbstaff-equipment\t	ransfer\DH_JM_Autosave	e_Test\3D_Time_TIF		
1		Data type :			TIF \$		
641	660	Compression :			Uncompressed 🗢		
	660	Shift to 16 bit : Bri	ightness optimized O Unprocessed		Å		
		Advanced image dat     Alternative data tvi	a configuration	)	*		
_							
						~	
							1 Million
		Counti	ng ∓) None ∓) ♥		and the state	20	1. 200

# INTRODUCTION

In this guide we describe the usuage of the Leica LASX auto-save functionality. By using auto-save, there is less risk of data loss during an extended image acquisition session and long file saving times after an extended imaging session can be avoided. Since auto-save does not generate standard Leica .lif files, we also show how the resulting files can be handled and visualized.

Here we show the workflow for multichannel, multiposition, 3D time series data.

#### Step 1 — Activate the auto-save function



- In "Open projects" at the bottom of the window you can find the autosave settings for the next project.
- Click on the "+" icon in the autosave settings to open the advanced configuration.
- Changes in the auto-save settings are only applied when generating a new project. So always make a new project when engaging or changing auto-saving.

			×				
1 f 1	<ul> <li>Project settings:</li> </ul>	0.*					
	Next project name :	3D_Time_Multi_TIF_Lossless					
	Auto-Save :	ON					
	Base data folder:	\\files.core.uzh.ch\zmbstaff-equipment\transfer\DH_JM_Autos Browse					
	Current working folder:	Browse					
	Current data path :	$\label{eq:linear} $$ $ $ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \$					
	Data type :	πε \$					
	Compression :	Lossless 🗧					
	Shift to 16 bit:  Brightness optimized  Unprocessed						
	<ul> <li>Advanced image dat</li> </ul>						
	<ul> <li>Alternative data typ</li> </ul>	e: LOF, Uncompressed, Shift to 16 bit: No					
	Data type :	LOF \$					

# Step 2 — Recommended auto-save settings

- We recommend the following settings for autosaving:
  - Data type: TIF
  - Compression: Lossless
  - Shift to 16 bit: Unprocessed
- Select a "Base data folder" with enough free space. The "Current working folder" can be left blank.
- The proprietary Leica .lof auto-save file format can be opened in LASX as well as in Imaris. However, the OME Bioformats reader can not read it. Due to this limited compatibility we recommend tif.

#### Step 3 — Perform your experiment and close the project



- Before starting your experiment make a new project with the updated auto-save settings.
- Perform your experiment. For multiposition experiments we recommend to use the LAS Navigator.
- Check that the data has been saved to your specified auto-save base data folder.
- If needed you can now still save your data in the .lif file format. However, this might take long.
- Close your project. Since all data has been saved, that should be very quick.
- Deleting the project will also delete the auto-saved data.

# Step 4 — Convert the resulting files in Imaris

Living Image	erviy Image I - Jak	De Imario Fale Converter				-					
Am		Enput Drag & Drop Piles, or click Button below to add Piles.	Duput			Select Files for Conversion. Co + + 2020_02_2  Organize + New folde	8,09,32,47~3D_Time_Multi_TIF_Lossless	1 → TileScan1 →		• 49 Search TieSea	12 P
Amira 2019. Anacond Navigato Arivis Arivis	2 Internet 930 Internet 4330	Address Soci preserver per reformer (social for a f	E (end/the:         2019/0004/2010/E (0.04.07 µundes).pp)         0           Data         Data         Data           Bester/End/terrenew.         Bester/End/terrenew.         Bester/End/terrenew.	Elected Input F	<b>t</b> oskopie		Name           Model/de           Partice 1           Partice 2           Partice 2           Partice 2           Partice 3           Parti	Date modified 28.02.2003 0.001 28.03.2003 0.001	Type File folder File folder File folder File folder File folder File folder File folder File folder File folder File folder Tiff image TIFF image	77 KB 48 KB 55 KB 17 KB	ů.
CellProfiler	LSX Mag View	Pedrarosa. Boot.		Canod Al		Without File na	Applient -00-201-00	28.0.2.202 09.33 28.0.2.202 09.33 28.0.2.202 09.33 28.0.2.202 09.33 38.0.3 30.00.00.23	TBF image TBF image TBF image TBF image	50.03 72.16 82.26 6. v 0 • All Files (1:5) Cipen	- Cancel

- Open Imaris File Converter.
- Choose an output folder.
- Select the directory where the tif files were saved.
  - When saving multi positions using the Navigator a **TileScan** folder should include all the tif files that were acquired.
  - Select the first file.

# Step 5 — Convert the resulting files in Imaris



- Press settings to check and adjust the correct dimensions.
  - In this case each position will be converted into an independent \*.ims file. (F split)
  - Each file contains a number of different time points (t),
  - a number of Z sections (Z),
  - and a number of channels.
- Start conversion.

(i) If you choose "Same folder as input file" copy the converted Imaris file to a new sub-folder.

#### Step 6 — Browse the resulting files in Imaris



- Open Imaris
- Add the folder containing your converted files in Arena by clicking "Observe Folder".
- Choose the directory where the converted files were stored.
- Here you will see a preview of your files for browsing your experiment.
- You can now open and explore each file individually.

# Step 7 — Open and browse the resulting files in FIJI



- You can also use FiJi to visualize and process your data.
- Open FiJi.
- Open files using File<Import<Bio-Formats.
- Open the folder containing the TileScan and select the first file.
- Select the options "Group files with similar names".

# Step 8 — Open and browse the resulting files in FIJI



- Confirm the dimensionalities. Eg.
  - Position 1--t00--Z00--C00.tif
  - Axis 1 10 Positions, Axis 2 10 time points, Axis 3 5 Z, Axis 4 4 channels
  - All the files will be loaded into a single hyperstack.

The different timepoints will need to be resorted.

 Make a new Substack with an increment to correctly display the time series. Eg. if you have 10 timepoints per position you will need to add an increment of 10.

#### Step 9 — Make new substack



- Image<Stacks<Tools<Make Substack</p>
- Make a new Substack with an increment to correctly display the time series. Eg. if you have 10 timepoints per position you will need to add an increment of 10.
- Image<Stacks<Tools<Make Substack</p>
- To display the 2 first positions:
  - 1-100-10,2-100-10