

ELYRA - Channel alignment

Correcting SIM images for chromatic shifts: This analysis step is only required for datasets containing more than one color channel.

Written By: z mbstaff

Step 1 — Channel alignment for SIM



- Collect data using SIM mode with the same imaging parameters as your sample.
- Acquire a z-Stack of TetraSpeckTM microspheres spread onto #1.5 coverslips and embedded in a hardening mounting media of your choice with a accurate step size in the relevant colors with SIM (identical settings as your experiment) and process the data with the 'SIM' processing tool.
 - (i) Alternatively, use the microspheres slide from Carl Zeiss Microscopy GmbH.
- This acquired stack will be used to calibrate the color offset of the system. Offsets will be stored in a template that later on can be used to correct the experimental data.
- Run the Structured Illumination processing as described in the guide.
- Calibrated color offsets are specific for the objective and filters. Therefore, a new calibration is required when objectives and filters are exchanged.

Step 2 — Obtaining the channel alignment transformation table from your bead data



- In the **Processing** tab select the Channel Alignment tool from the list of Methods.
- In the Method Parameters tab select the file that you obtained from the Structured Illumination processing of your bead stack by opening the file in ZEN and clicking Select. The filename of the selected dataset will appear next to the button.
- In the Channel Alignment options make sure Fit is checked and select Affine, Lateral or Parabolic from the dropdown menu.
 - Lateral is used to correct for linear lateral and axial drifts, that is in all directions. Linear alignments are performed in the x, y and z direction.
 - Affine is used to correct for linear lateral (X, Y) and axial (Z) drifts, rotational offsets in the image plane in x and y (RX, RY) as well as a 1st order stretching in the image plane in x and y (SX, SY).
- Press **Apply** near the top of the screen to start the processing (takes some time to run).
- In the resulting image all colour channels should be properly aligned now. Save the transformation table on the left as a ".bin" file by clicking the Save-button.

Step 3 — Applying the channel alignment to your SIM images



- In the Processing tab select the Channel Alignment tool from the list of Methods.
- In the Method Parameters tab select the Structured Illumination output image of your sample that you want to correct for chromatic aberrations by opening it in ZEN and clicking Select. The filename of the selected dataset will appear next to the button.

Step 4 — Applying the channel alignment to your SIM images

Method Para	meters test_Structured Illumina	ation	Select	
Output Image	test_Structured IIIu	stion_Channel A	put 2	
ID Laser 0 405 1 488 2 561 3 5642	X Y 0 0 -0.345i 0.2543 -0.314i 6.3288 -0.996i 0.9585	RX RY 0 0 8.1008 5.9713 -0.000 -0.000 0.0003 -1.437	SX 0 1.6792 -0.000 0.0006	
CH La 17V2-73 48 17V2-72 56 Set Defaults	ser ID 8 1 • 1 2 • Get Defaults	Save Lo	bd	

- Unselect the 'Fit' check box
- Then click on the Load-button under the empty transformation table, select the bin file you saved earlier and press Open. Now the table should display the values for the colour shifts between channels.
- .bin file is used to compensate for any chromatic aberrations and offsets between channels in SIM.
- In the table below make sure the wavelength channel IDs correspond to the wavelengths of the channels in the transformation table above.

- Press Apply to apply the transformation table to your sample image.
- Save the resulting file in .czi format.
 ZEN will automatically append
 "Channel Alignment" to the output filename.
- Refractive index mismatches don't affect the correction template as the beads are on the coverslip. If the sample is fixed to the coverslip, the correction is reliable regardless of embedding differences.
 However, if fixed on the slide, mismatches may impact accuracy, requiring beads to be embedded in the same medium as the sample.