


# Spinning Disk - Visitron (Irchel): Shut down

Written By: Jana Döhner




## Step 1 — Cleaning and closing software

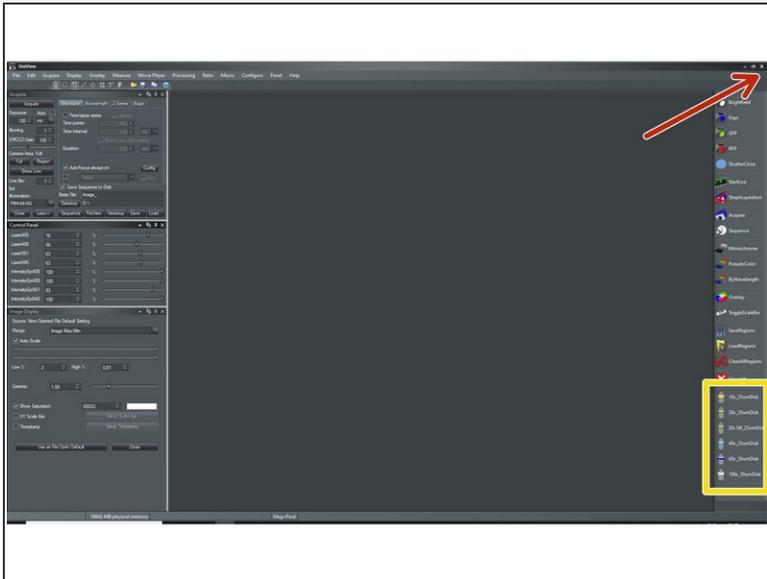


- **Clean** the used immersion objectives using the available **KIMTECH wipes** and **100% EtOH**.
- **Close** the software by crossing it out

## Step 2 — Stage Insert

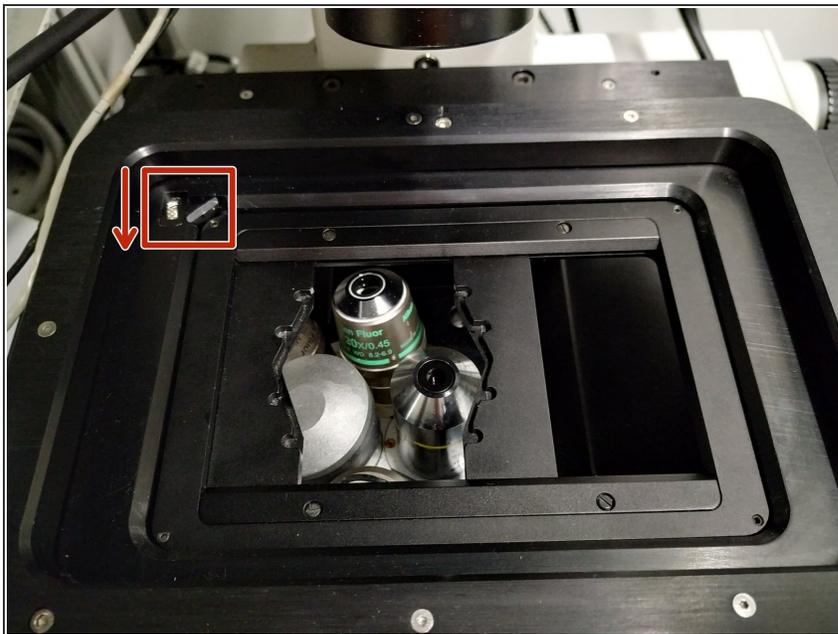

- **Remove** the used **stage insert** by releasing the screw (arrow direction indicates how to loosen the screw) and **stow away** accordingly

### Step 3 — Log-Off or Shut Down

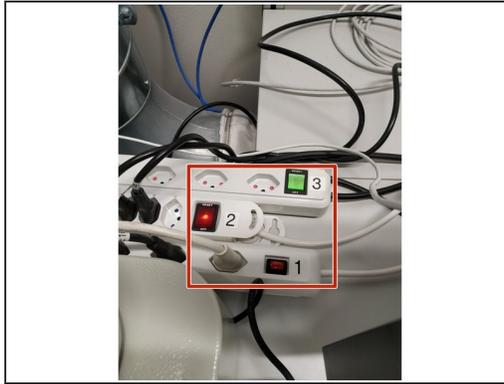

- i** Check in the **reservation system** if there is another booking shortly after you.
- If **"YES"** keep the laser keys ON, and just sign-out from the PC.
  - If **"NO"** shut-down the PC and turn OFF all laser keys (4).
  - After the PC has been shut-down switch OFF all power supplies (3, 2 and 1)
  - **Cover** the microscope.